

Selection of Stable Transfected Cell Lines

Background:

Once the cells have been successfully transfected, the next step is to seed and select the transfected cell line in a single 96-well plate to select pure colonies by limited dilution as outlined below:

Protocol:

- 1. Seed the transfected cells in 96-well plates in 10% conditioned medium:
 - a. Dilution conditions:
 - 2x96 well plate with 0.1 cell per well
 - 2x96 well plate with 0.5 cell per well
 - 2x96 well plate with 1 cell per well
- 2. Incubate the cells for 24h.
- 3. Remove medium and add conditioned selection medium containing selection antibiotic at the predetermined concentration required for your cell line. Incubate 96-well plates at 37°C with CO2.
- 4. Check the plates every day for colonies. Colony formation depends on proliferation rate of the cell line and can take anywhere from 3 days to 1 week.
- 5. Refresh selective medium every 3-4 days until colonies appear.
- 6. Select the wells with only one single colony. Make sure colonies are not growing in clumps as they will be less sensitive to the antibiotic.
- 7. When a well contains a single colony, transfer the colony to a 24-well in selection medium and so on until you have enough cells for freezing and storage in liquid nitrogen. Use the appropriate antibiotic concentration as determined from the kill curve.

QC

Seed 24-wells with insert for an immunostaining to determine percentage of cells expressing the gene of interest to be able to identify a 100% pure clone. You can also use Western blotting, flow cytometry or another technique depending on the cell line used.

Seed 24-wells with insert and determine the expression level of the gene of interest by immunostaining.

- 1. Grow cells on insert in a 24-well plate until well has confluent growth.
- 2. Remove medium and wash cells with 1X PBS.



- 3. Fix cell with methanol or paraformaldehyde and wash with 1X PBS.
- 4. Add primary antibody in 24-well against protein of interest and incubate at 37°C for 1 hour (depending on antibody).
- 5. Wash cells with 1X PBS.
- 6. Add secondary antibody in 24-well plate and incubate at 37°C for 1 hour (time depends on antibody type).
- 7. Wash cells with 1X PBS.
- 8. Remove insert from 24-well plate and affix to microscopy slide with nail polish or other appropriate adhesive.
- 9. Determine the percentage of transfected cells with fluorescence microscope.