



Restriction Enzyme Afl II



Cat.#	Size	Conc.
FG-AflII	2,000 units	4 units/μl

Store at -20°C

Supplied with: 10X FastGene® Buffer IV (FG-REB4)
10X FastGene® FastCut Buffer (FG-REBHF)
6X DNA Loading Buffer
Sterile water

Recognition site



For Research Use Only. Not for use in diagnostic procedures.

ISO9001

Source: *Anabaena flos-aquae*

Reaction conditions

1X FastGene® IV 37°C
1X FastGene® FastCut Buffer, 37°C

FastGene® FastCut Buffer

FastGene® restriction enzyme can cut substrate DNA in 5-15 with FastGene® FastCut Buffer.

1X FastGene® Buffer IV

20 mM Tris-acetate (pH 7.9 at 25°C)
50 mM potassium acetate
10 mM magnesium acetate
100 μg/ml BSA

Unit definition

One unit is defined as the amount of enzyme required for complete digestion of 1 μg bacteriophage λ (Hind III digestion) at 37°C for 1 hr in 50 μl reaction mixtures.

Quality control

- Unit definition assay
- Overdigestion assay
- Endonuclease assay
- Extreme pure assay

Dilution buffer:

FastGene® Diluent A

Heat Inactivation

Afl II can be inactivated at 65°C for 20 min.

Methylation sensitivity

dam methylation: Not sensitive
dcm methylation: Not sensitive
CpG methylation: Not sensitive

Prolonged incubation

A minimum amount of enzyme required to digest 1 μg substrate DNA for 16 hr; 0.13 U.

Relative activity in FastGene® Buffers

FastGene® Buffer I: 75%
FastGene® Buffer II: 100%
FastGene® Buffer III: 75%
FastGene® Buffer IV: 100%
FastGene® FastCut Buffer: 100%

Note

It is not sensitive to *dam*, *dcm* or CpG methylation. Less than 50% of cleaved fragments are ligated in 20 μl reaction mixtures containing 100-500 units of T4 DNA ligase.

Standard reaction condition

- Normal protocol

Component	Final Conc.	Volume
Substrate DNA	1 µg	X µl
10X FastGene® Buffer IV	1 X	5 µl
Afl II	4 unit	1 µl
Sterile water	up to 50 µl	
→ Incubate at 37°C for 1 hr		

- Fast protocol

Component	Final Conc.	Volume
Substrate DNA	1 µg	X µl
10X FastGene® FastCut Buffer	1 X	5 µl
Afl II	4 unit	1 µl
Sterile water		up to 50 µl
→ Incubate at 37°C for 15 min		

※ We recommend 5-10 units of enzyme per μg DNA and 10-20 units for genomic DNA in a 1 h digest



Restriction Enzyme Afl II



Cat.#	Size	Conc.
FG-AflII	2,000 units	4 units/μl

Store at -20°C

Supplied with: 10X FastGene® Buffer IV (FG-REB4)
10X FastGene® FastCut Buffer (FG-REBHF)
6X DNA Loading Buffer
Sterile water

Recognition site



For Research Use Only. Not for use in diagnostic procedures.

ISO9001

Source: *Anabaena flos-aquae*

Reaction conditions

1X FastGene® Buffer IV 37°C
1X FastGene® FastCut Buffer, 37°C

FastGene® FastCut Buffer

FastGene® restriction enzyme can cut substrate DNA in 5-15 with FastGene® FastCut Buffer.

1X FastGene® Buffer IV

20 mM Tris-acetate (pH 7.9 at 25°C)
50 mM potassium acetate
10 mM magnesium acetate
100 μg/ml BSA

Unit definition

One unit is defined as the amount of enzyme required for complete digestion of 1 μg bacteriophage λ (Hind III digestion) at 37°C for 1 hr in 50 μl reaction mixtures.

Quality control

- Unit definition assay
- Overdigestion assay
- Endonuclease assay
- Extreme pure assay

Dilution buffer:

FastGene® Diluent A

Heat Inactivation

Afl II can be inactivated at 65°C for 20 min.

Methylation sensitivity

dam methylation: Not sensitive
dcm methylation: Not sensitive
CpG methylation: Not sensitive

Prolonged incubation

A minimum amount of enzyme required to digest 1 μg substrate DNA for 16 hr; 0.13 U.

Relative activity in FastGene® Buffers

FastGene® Buffer I: 75%
FastGene® Buffer II: 100%
FastGene® Buffer III: 75%
FastGene® Buffer IV: 100%
FastGene® FastCut Buffer: 100%

Note

It is not sensitive to *dam*, *dcm* or CpG methylation. Less than 50% of cleaved fragments are ligated in 20 μl reaction mixtures containing 100-500 units of T4 DNA ligase.

Standard reaction condition

- Normal protocol

Component	Final Conc.	Volume
Substrate DNA	1 µg	X µl
10X FastGene® Buffer IV	1 X	5 µl
Afl II	4 unit	1 µl
Sterile water	up to 50 µl	
→ Incubate at 37°C for 1 hr		

- Fast protocol

Component	Final Conc.	Volume
Substrate DNA	1 µg	X µl
10X FastGene® FastCut Buffer	1 X	5 µl
Afl II	4 unit	1 µl
Sterile water	up to 50 µl	
→ Incubate at 37°C for 15 min		

※ We recommend 5-10 units of enzyme per μg DNA and 10-20 units for genomic DNA in a 1 h digest